


Lambda *EcoRI* Ladder

Catalogue Number: 1LLE-500
Volume: 5 x 1 ml for 500 loadings

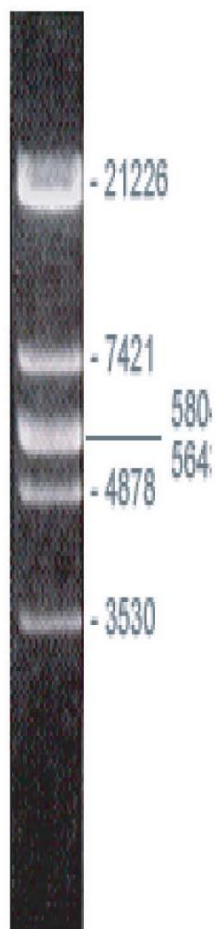

The λ *EcoRI* DNA Ladder is prepared by restriction digestion of Lambda DNA to completion with *EcoRI*, followed by heat inactivation of the enzyme.

The resulting 6 fragments range from 3.53 kilobases to 21.23 kilobases in size. (3530, 4878, 5643, 5804, 7421, 21226 bp)

The λ *EcoRI* DNA Ladder is supplied ready to use in loading buffer containing MZ Blue Dye.

Load 10 to 15 μ l onto the gel (340 to 500 ng)

Recommended concentration of agarose 0.7 to 1.7%.