


Microzone Ltd

Lambda *Hind*III Ladder

Catalogue Number: 1LLH-500

Volume: 5 x 1 ml for 500 loadings

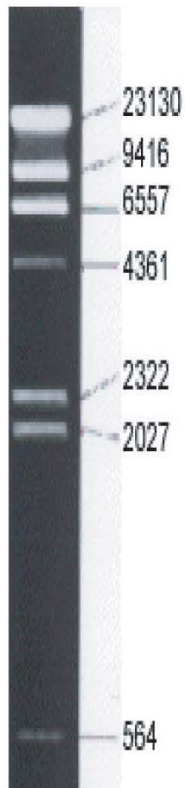

The λ *Hind*III DNA Ladder is prepared by restriction digestion of Lambda DNA to completion with *Hind*III, followed by heat inactivation of the enzyme.

The resulting 7 fragments range from 564 base pairs to 23.13 kilobases in size. (564, 2027, 2322, 4361, 6557, 9416, 23130 bp)

The λ *Hind* DNA Ladder is supplied ready to use in loading buffer containing MZ Blue Dye.

Load 10 to 15 μ l onto the gel (340 to 500 ng)

Recommended concentration of agarose 0.7 to 1.7%.



Store at +4 or -20°C

For Research Only