

BetterBuffer Sequence Enhance/Dilution Buffer

Specially formulated to be used with **ABI BigDye® Mix (versions 1.0 and 3.0)**

Protocol

- Dilute the ABI BigDye Mix with BetterBuffer to achieve the desired concentration*
- Mix by pipetting
- Sequence according to protocol
- Remove excess dye

*1 in 6, 1 in 12, 1 in 15 or higher. The choice of dilution is dependent on size, quantity and quality of the DNA template. Please contact Microzone for answers to any questions
BigDye is a registered trademark of Applied Biosystems

Related Products

microCLEAN = The easy way to clean-up your PCR fragments

Just Water = Molecular grade water in handy sizes

Supplied as a clear solution in multiple of **0.3, 1.5, 25 & 250 ml aliquots**
Store at **-20°C**

For Research Only